# LAVAJOS DE EMERGENCIA EMERGENCY EYE WASHERS

## Nahita

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
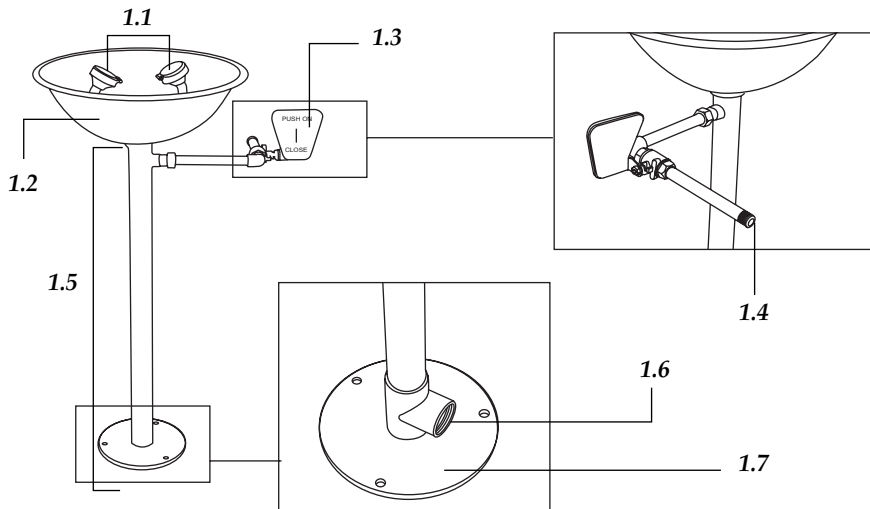
CODE.:64000759

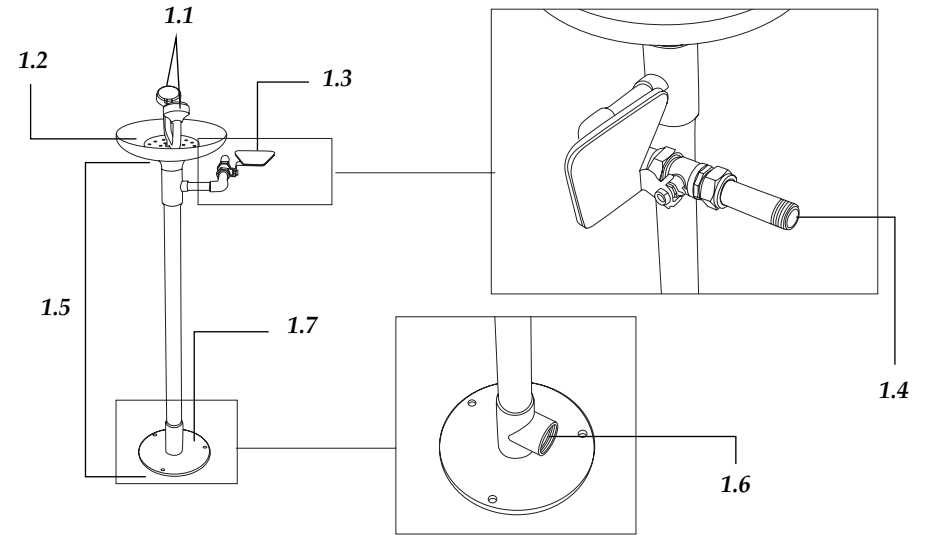**4. INSTRUCTIONS OF USE**

*Note: Eye washer must be use only in case of emergency, not for a daily use. Activate the eye washer at least weekly to assure that it will operate properly when needed.*

- ◆ To activate the sprays push the lever clockwise.
- ◆ Keep eyes opened with the help of hands so as the damaged zone is abundantly washed with water.
- ◆ To stop water flow, press the lever counter-clockwise.

*Note: The time of wash must be adequate to the type of damage produced. The emergency eye washers can reduce the degree of damage produced but they are never substitutive of a medical treatment. Apply the appropriate medical treatment.*

We will only accept the product return within 15 days after delivery provided it comes in its original wrapping.

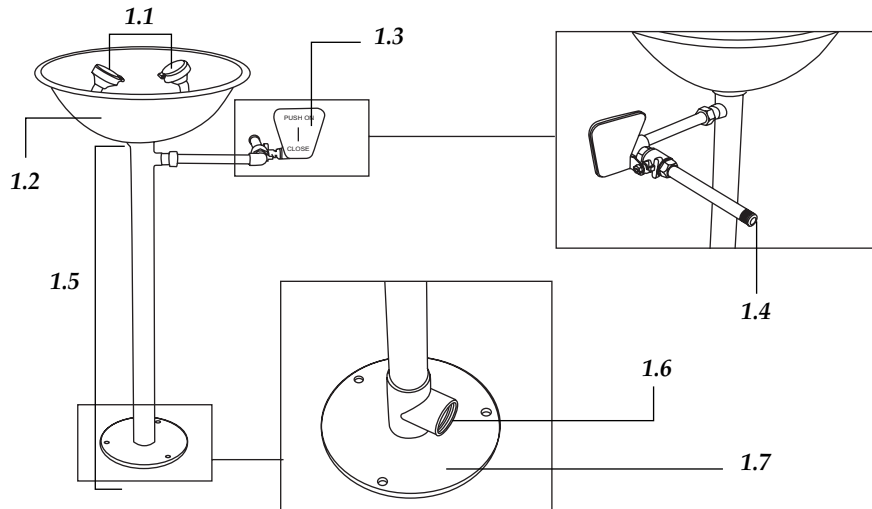
### Installation

**IMPORTANT:** The place chosen to install the eye washer must be adequate, so as it is perfectly visible and accessible from the region where the risk of accident exists. The distance must not be higher than 15 m o 10 s walking from the zone of risk.


- ◆ Wrap all threads with teflon tape or similar to assure all joints are perfectly sealed and avoid water leaks. Then, assemble the different pieces of the eye washer according to the diagram shown:
- ◆ Screw the main body in the base of the eye washer so as water outlet is downwards.
- ◆ Install the tank and then the two sprays head.
- ◆ Put the starter lever so as it opens the valve when being pushed clockwise.
- ◆ Connect the eye washer to the water inlet and outlet.

- |                   |                  |
|-------------------|------------------|
| 1.1 Sprays        | 1.5 Main body    |
| 1.2 Tank          | 1.6 Water outlet |
| 1.3 Starter lever | 1.7 Base         |
| 1.4 Water inlet   |                  |

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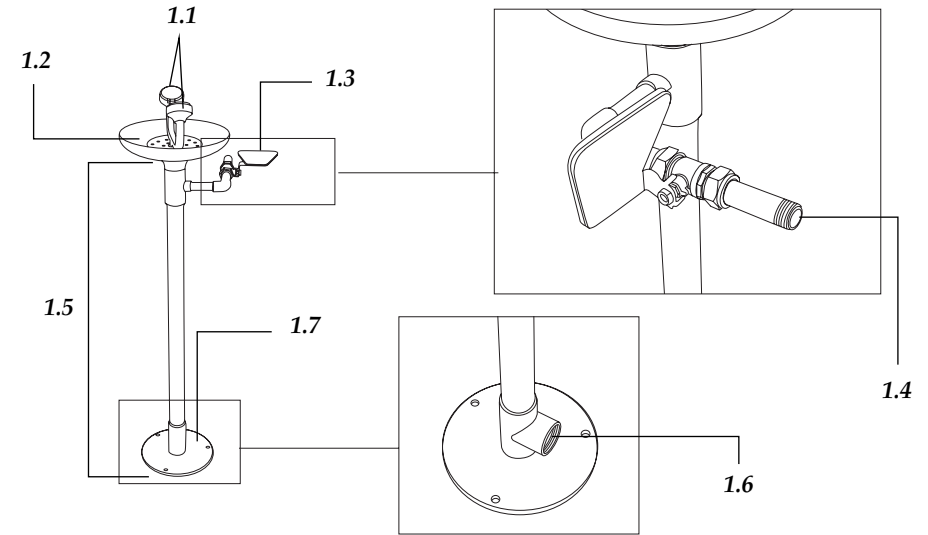
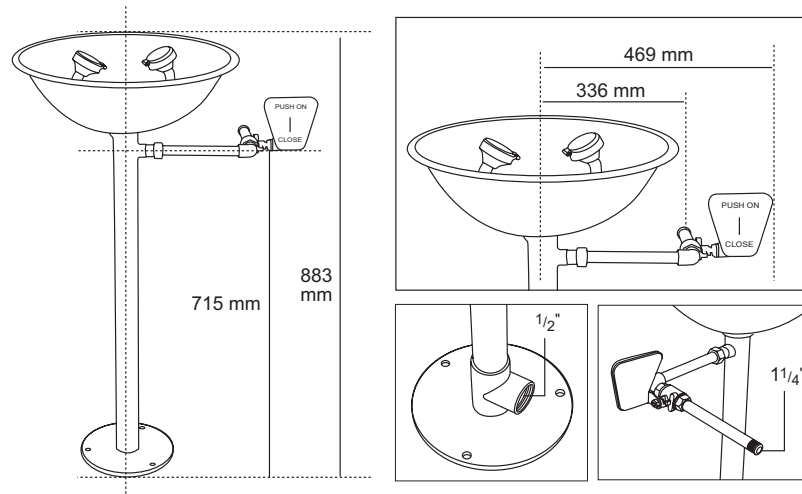
REF.:64000759

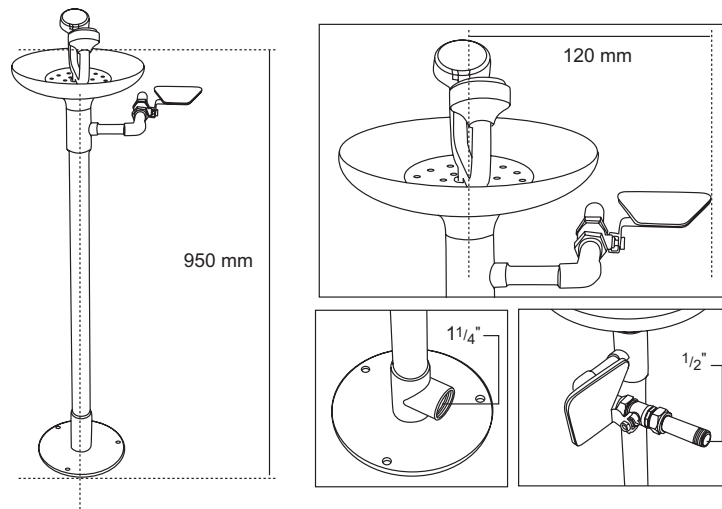**1. USES**

Emergency showers and eye washers are a necessary security measure with high utility in most of companies and laboratories. They can reduce the damage caused by the splashing in the eyes, face and the body with acids, bases, organic compounds and other toxic or corrosive components, as well as in case of fire. It is the easiest and cheapest way of decontamination in case of accident.

**2. TECHNICAL SPECIFICATIONS**

Pressure			Flow (L/min)	Temperature
Nominal pressure	Seal pressure	Working pressure	≥ 11.4	Tepid water
0.4	0.45	0.20-0.40		

- ◆ Made of high quality materials: stainless steel (model 64000359) or anticorrosive ABS plastic tank and PVC-coated steel pipe (model 64000759).
- ◆ The base can be fixed to the floor to assure a high stability.
- ◆ High flow sprays of low pressure water to avoid damages. Protected from dust and other contaminants with caps that are removable without the intervention of user.
- ◆ The unit is designed with enough space to keep eyes opened with the help of hands while water flows towards the eyes.
- ◆ The valve is automatically opened when pressing the starter lever; it allows a permanent water flow until it is closed manually by pressing the lever to the opposite direction.

**3. INSTALLATION**

Preliminary inspection

- ◆ Unwrap the eye washer, take off all the protective items and make sure that it does not present any damage because of the shipment. In case the eye washer presents any damage tell it immediately to your transport agent or dealer so that they can make the claims in the correct time limit.
- ◆ Please keep the original wrapping; you will always need it for returns enclosed with all the accessories supplied.
- ◆ Please check that all the parts are enclosed:
  - Head with two sprays
  - Tank
  - Main body
  - Base
  - Starter lever
  - User's manual

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958 412 886

629 636 705

<http://www.cromakit.es/>

**Calle Tucumán 8 Nave B 18200 Maracena (Granada)**